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MECHANIC

SHOW THE STYLE OF MASTER

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Contents

1. Before use	2
1. 注意事項.....	2
2. Nomenclature	3
2. 各部件名称	3
3. Operation	4
3. 各部件安装	4
4. Configuration	8
4. 配置表	8
5. Technical parameter	9
5. 技术参数.....	9
6. Troubleshooting	10
6. 故障处理.....	10

1 Before use 注意事项

1-1 NOTICE

- 1) Microscope ought to be placed in a dry and clean place. Do not expose the microscope in the sun directly. Avoid high temperature and violent vibration.
- 2) As microscope is a precision instrument, handle with care, avoiding impact or abrupt movement during transportation.
- 3) To keep the image clear, do not leave fingerprints or stains on the surfaces of the lens.
- 4) Never turn the left and right focusing knob in the adverse direction at the same time, otherwise the microscope will be damaged.
- 5) Hold the camera with one hand for fearing of falling when you take the films out of the big camera.

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1-2 MAINTENANCE

- 1) All lenses must be kept clean. Fine dust on the surface of the lens should be blown off with hand blower or wiped off gently with a soft lens tissue; Fingerprints or oil marked on it should be wiped off with a tissue moistened with a small amount of xylene or a 3:7 mixture of alcohol and ether.
- 2) Never use the organic solution to clean the other surface (especially the plastic surfaces). If necessary, please choose the neutral detergent.
- 3) Do not take the microscope apart for fearing that it is damaged.
- 4) After using, cover the microscope with the dust-cover provided and store it in a dry and clean place free from moisture to prevent rust.
- 5) To keep the performance of the microscope, please check it periodically. The detail can be gotten from the agent nearby.

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
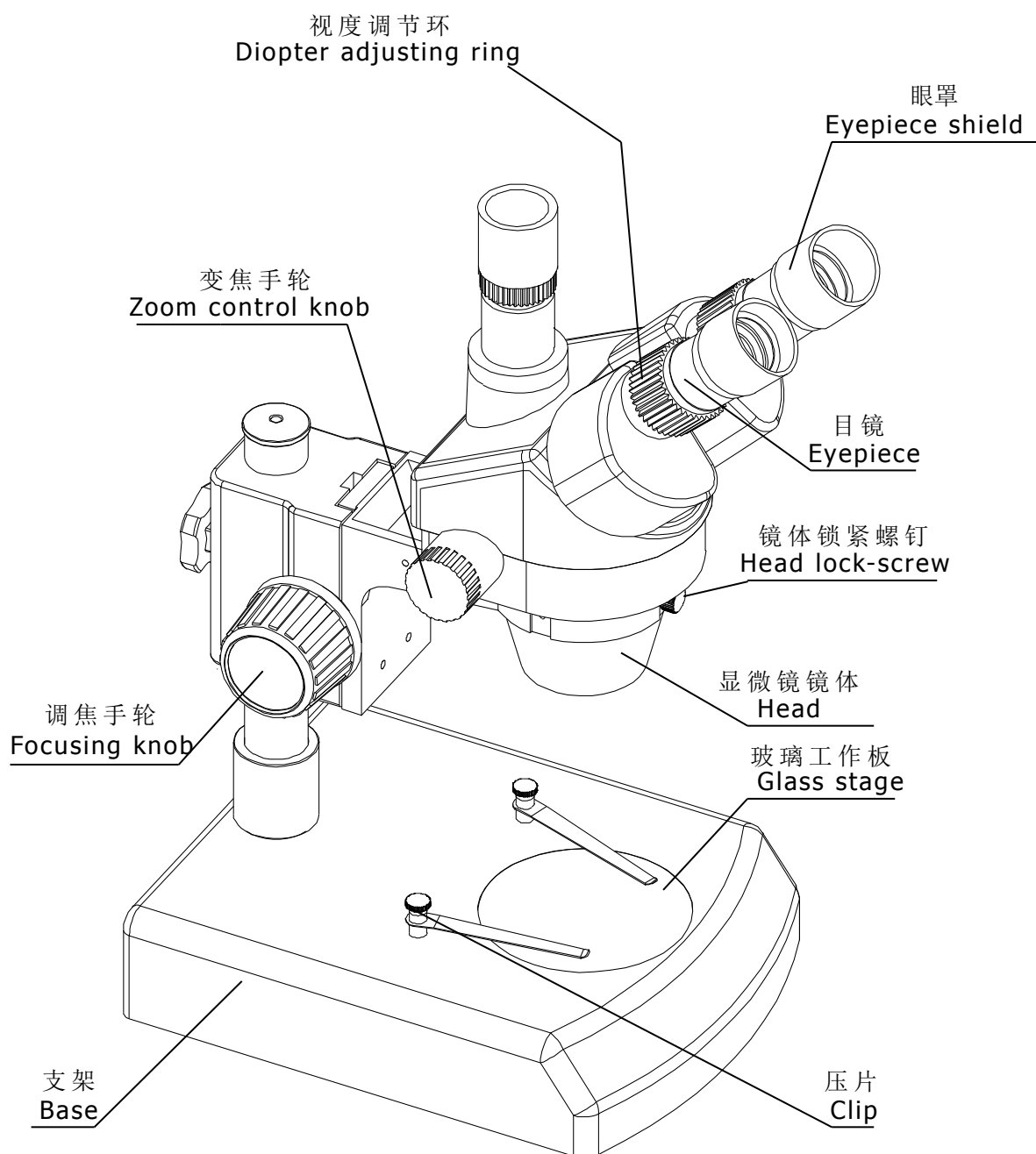
3 Operation 操作

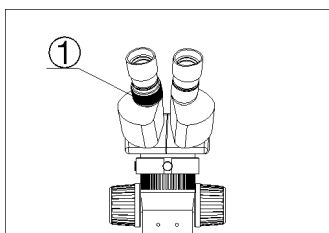3-1 Place the specimen

Place the specimen in the middle of the glass workbench. If it is necessary to hold down the specimen, choose a suitable lighting method to illuminate the specimen.

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3-2 Viewing adjustment and focus

Observe through the eyepiece on the right, rotate the focusing handwheel to make the specimen image clear observe through the left eyepiece, rotate the left diopter adjustment ring to see the specimen clearly

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3-3 Adjust interpupillary distance

Hold the left and right prism boxes with your hands, and rotate the left and right prism boxes in the direction shown by the arrow until you feel comfortable in binocular observation.

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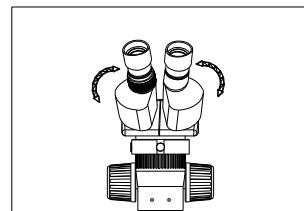
3-4 Use of eye mask

For people who do not wear glasses, hold the diopter adjustment ring with your hands so that it does not rotate. By rotating the eyepiece, the eyecup can fit the observer's eye better.

For people with glasses, you can take off the goggles and observe

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3-5 Installing and removing the reticle

Unscrew the reticle holder from the eyepiece. Clean the reticle, place the engraved side of the reticle down, and install it into the reticle holder. Screw the reticle holder with the reticle into the eyepiece holder until it is tightened. To disassemble the reticle, unscrew the reticle holder from the eyepiece, take out the reticle, wrap it in clean paper, and store it

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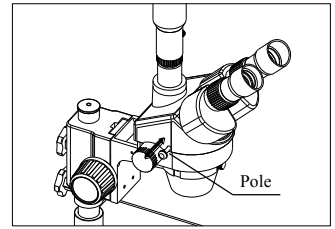
3-6 choose the optical system

You can alternate the binocular observation and video capture by pushing or pulling the pole. You can attain binocular observation by pushing the pole inside, or attain video capture by pulling it outside. No matter what optical system is chosen, push or pull the pole thoroughly.

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
3-7 Mount the photo eyepiece and the PK-mount adapter

Put the photo eyepieces socket of the trinocular. Connect the PK-mount adapter with the photo eyepiece, and then tighten the lock-screw.

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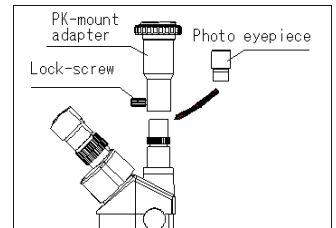3-8 Adjust the CTV

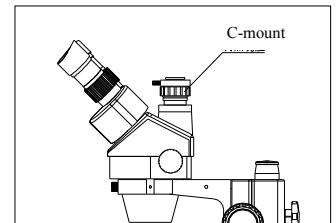
Adjust the CTV to a suitable position by rotating C-mount.

Note: The range of the adjustment: 1~2mm in general.

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3-9 Connect the Digital head With the Monitor or TV set

1) Plug one end of the PVA cable into the socket of the digital head.

2) Plug the C-VIDEO or S-VIDEO of the PVA cable into the correct socket of Monitor (TV set).

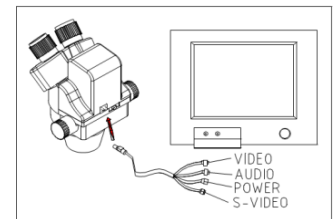
3) Connect the 12V DC power with the power socket of the PVA cable.

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3-10 Appear the image on the Monitor or TV

1) Connect the power supply and then turn on the Monitor or TV.

2) For the monitor, the connect sign model must be chosen (C-video or S-video) and for TV, the channel must be set to the video channel.

3) Pull pole out and adjust the focusing knob and then the image will appear on the screen clear

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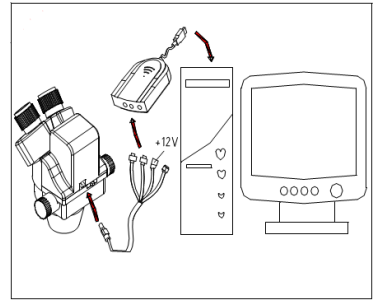
3-11 Connect with the computer

- 1) Plug one end of the PVA cable into the socket of the digital head.
- 2) Plug one of the C-VIDEO or S-VIDEO into the A/D board.
- 3) Plug the USB of the A/D board into the USB socket of the computer.
- 4) If your computer is mounted capture card, you can connect the C-VIDEO or S-VIDEO with the computer directly.
- 5) Connect 12V DC power with the power socket of the PVA cable.

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3-12 Appear the image on the computer

- 1) Turn on the power supply and let the computer work.
- 2) Install the software and the driver of the A/D board.
- 3) Double click the icon of the software, and then the video window will appear. You can set the size of the window according to your linking
- 4) Draw out the pole and adjust the focusing knob, and then the image will appear on the computer screen clearly.
- 5) If no image or the image without color, it may be because the model of the input signal does not match the output signal of CCD or the model of C-VIDEO/S-VIDEO is not correct. The detail of operation refers to «Software operation manual»

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3-13 Appear the image on the computer and the Monitor synchronously

- 1) Do step 4-12 and step 4-14 to connect the computer and the Monitor.
- 2) Operate step 4-13 and step 4-15, we can make the image appear on the computer and Monitor at the same time.

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3-14 Adjust the image

- 1) Put the base, stand and digital head correctly, then fix the lock-screw tightly.
- 2) Put the object on the base stage.
- 3) Observe the object through the eyepiece and adjust the focusing knob to make the image of the object clearly.
- 4) Move the digital head or the object gently to adjust the image agreeing with observer.

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3-14 Brief instruction for the software

- 1) The program design of the software is up to date, and the Chinese/English interface can berth powerful delineation bar which be used much conveniently and rapidly. You can finish most of analyze work only to click the mouse.
- 2) Can afford many powerful area choosing tools which can analyse any area your linking at will, such as adjusting hue and image, dealing with mathematical morphology, image matching, texture analyses, character identify and so on.
- 3) Geometry character measuring function, automatically analyzing function such as slightness body, grain body, line body and so on. The outcome can be kept in data and can be made into chart and so forth.

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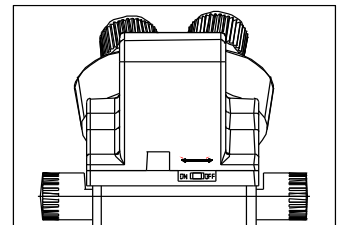
3-15 Use the white balance

- 1) The CCD has auto white balance when the white balance switch is on "ON".
- 2) Please put the switch on "ON" in general. Let the switch be "OFF" only in special, for example, observing the red cell, otherwise the color of red cell will be adjusted into white.
- 3) If you want to observe another single color, please let the switch be "ON" again when you finish the observation, and put the switch on "OFF" again after auto balance, or the color of the image will be distortion.

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3-16 Use of Base work board

You can lift the base working plate directly from the recess with your fingertips.

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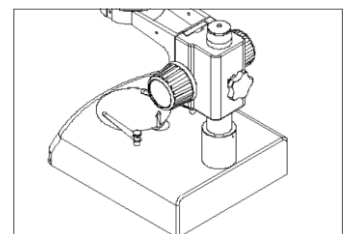
3-17 Tuning tightness

To adjust the tightness, you can hold one of the wheels with your hand and rotate it. The degree of tightness depends on the direction of rotation of the wheel: clockwise tight, counterclockwise loose. The tightness is suitable to prevent the microscope body from sliding down during the observation process, and it also makes the hand feel comfortable

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4 configuration 配置表

4-1 MC75 series configuration

Note: The items marked "O" included and others for option

配置表格	Specification 规格	75T-B1	75T-B3	75T-iMatX	75T-iSlide	75T-iMat mini	75T-iMatPRO	75T-L1	75T-L2	75T-L3	
	pics 图片										
Eyepieces 目镜	10X20	O	O	O	O	O	O	O	O	O	
	10X22										
	10X23										
	20X10										
Tri-ocular 三目	75T	O	O	O	O	O	O	O	O	O	
	MCN-0.3X										
	MCN-0.5X										
	MCN-0.7X										
Auxiliary objective 辅助物镜	MCN-2X										
	focusing arm 托架	A1	O	O	O	O	O	O	O	O	O
		B1	O								
		B3		O							
Base 底座		iMatX			O						
	iSlide				O						
	iMat mini					O					
	iMatPRO						O				
	L1							O			
	L2								O		
	L3									O	
	L3									O	

MCN-0.3X

MCN-0.5X

MCN-0.7X

MCN-2X

B1

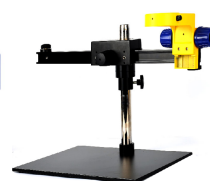
B2

L1

L2

L3

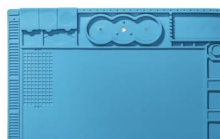
iMat mini

iMatPRO

iMatX

V56

iSlide

5 Technical parameter

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5-1 MC75T

MC75T Standard configuration 标准配置			Auxiliary objectives 辅助物镜					
			0.5X		1.5X		2X	
Eyepiece 目镜	Working distance 工作距离 100mm		Working distance 工作距离 165mm		Working distance 工作距离 45mm		Working distance 工作距离 30mm	
Type 型号	Magnification 放大倍率	Field of view 视场范围	Magnification 放大倍率	Field of view 视场范围	Magnification 放大倍率	Field of view 视场范围	Magnification 放大倍率	Field of view 视场范围
20X/10	14X	14.3	7X	28.6	21X	9.5	28X	7.1
	90X	2.2	45X	4.4	135X	1.5	180X	1.1
10X/20	7X	28.6	3.5X	57.2	10.5X	19	14X	14.2
	45X	4.4	22.5X	8.8	67.5X	3	90X	2.2
10X/22	7X	31.46	3.5X	62.92	10.5X	20.9	14X	15.62
	45X	4.84	22.5X	9.68	67.5X	3.3	90X	2.42
10X/23	7X	32.89	3.5X	10.12	10.5X	21.85	14X	16.33
	45X	5.06	22.5X	65.78	67.5X	3.45	90X	2.53

★Working distance is fixed regardless of the magnification factor.

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★Total mag. = Zoom mag. X Eyepiece mag. X Auxiliary objective mag.

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$$\text{Diameter of field of view (mm)} = \frac{\text{Field number of eyepiece 目镜视场}}{\text{Zoom mag. X Auxiliary objective mag. 物镜放大率 X 辅助物镜倍率}}$$

★Photo adaptor mag. = Zoom mag. (X Auxiliary objective mag.) X Eyepiece mag.

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6 Troubleshooting

The performance of the microscope can't be made fully because of unfamiliar using, this table will give some advices.

Trouble	Cause	Remedy
1、 Double images	Interpupillary distance is not correct	Readjust it
	Diopter adjustment is not correct	Readjust it
2、 Dirt appears in the field of view	Magnification of each eyepiece is not the same size	Mount the same size eyepiece
	Dirt on the specimen	Clean the specimen
3、 Image is not clear	Dirt on the surfaces of eyepiece	Clean the surface
	Dirt on the surfaces of the objectives	Clean the objectives
4、 Image is not clear while the focus changing	Diopter adjustment is not correct	Readjust the diopter
	Focus is not correct	Readjust the focus
5、 The focusing knob is not smooth	The focusing knob is too tight	Loosen it to a suitable position
6、 The image is obscure because of the head slipping down by itself during observation	The focusing knob is too loose	Tighten it to a suitable position
7、 Incision image appears in the field of view or of video view	The pole is not in correct position	Pull or push it to the correct position
8、 The image on the monitor is not clear when the focusing knob is turned.	The focus of video is not correct	Readjust the focus of video to a correct position
9、 Eyes fell tired easily	Diopter adjustment is not correct	Adjust the diopter
	Brightness of light is not correct	Adjust the brightness
10、 Bulb does not work when the switch is on	No power supply	Check the connection with the power supply
	The bulb was not inserted correctly	Insert it correctly
	Bulb is wrong	Replace with a new one
11、 Bulb is burned out suddenly	Use the wrong bulb	Replace with a correct one
	The voltage is too high	Control the voltage Eg: use voltage regulator
12、 Brightness is not enough	Use a wrong bulb	Replace with a correct one
	The voltage is too low	Increase the input voltage
13、 The bulb flickers or the brightness is unstable	The bulb will burn out soon	Replace with a new one
	The bulb was not inserted correctly	Insert it correctly

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